

# Pads / Plates cleaning

12/2012

- 45 min in Bleach.
- Rinse 3x with desalted water.
- 15 min. in 70% EtOH.
- Short in 100% EtOH.
- For Plates: bring the plates under the hood (transport-box) for drying.
- For Pads: put the pads in the boxes (laterally) with the paper already in and let them dry in the hood with the UV-light on.

Solution:

**Bleach:** dilute the bleach from 12% to 10% in the big container (10L volume is marked with a line).

Reference: Carl Roth (Natriumhypochloritlösung 12% Cl)  
Art Nr-9062.1.